

10 September, 2005

Instructions for JEOL MS-MP9020D data system:

- 1) On the JEOL workstation, from the Data manager or Data drawer window selected the raw file data.
- 2) Open the GCMS raw file containing the mass spectrum you wish to submit to MaSC.
- 3) Display the mass spectrum of interest and select Export to file - JEOL-DX (Right hand mouse button from the bar on top of the spectrum, Export to file - JEOL-DX). The range displayed on the window will be the range exported. The exported file will have a .jsp extension Retrieve the spectrum.
- 4) Open the terminal window (Dtterm). Go to the JEOL-DX export directory:
  - a. **cd /ms7801/Export/JEOL-DX/xms (setup dependent)**
- 5) Copy the .jsp file to the ftp directory: **Cp "yourfilename" /home/ftp/pub (setup dependent)**
- 6) Import the file into a PC using the ftp.